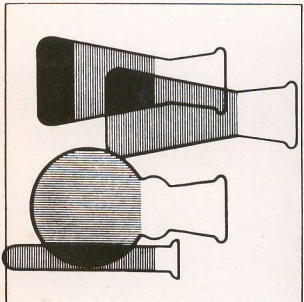
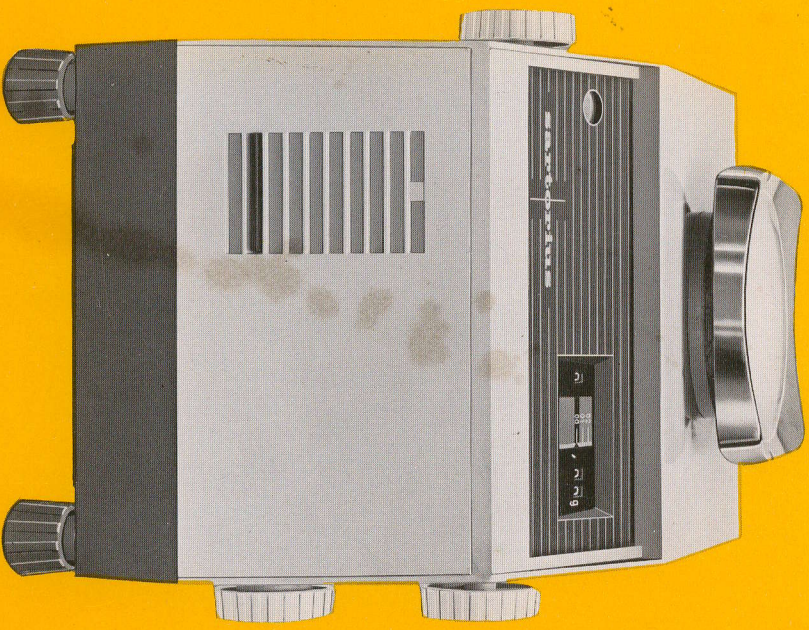
Precision Balances  
Balances de Précision  
Balanzas de Precisión  
Präzisionswaagen


# 2250

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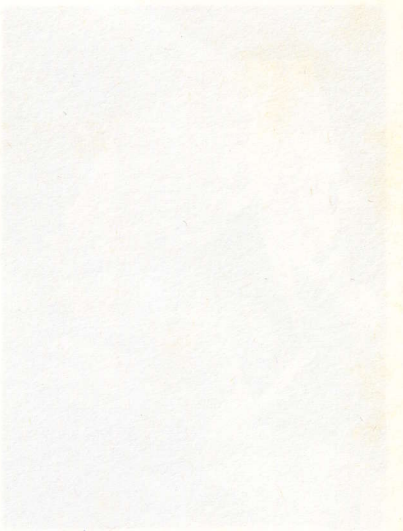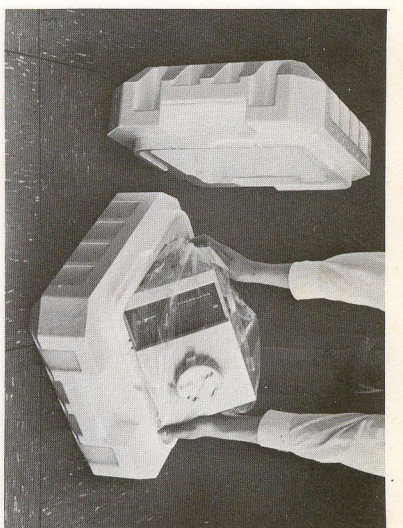
Remove Styrofoam insert from carton.

2

Open Styrofoam box and lift balance out.

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

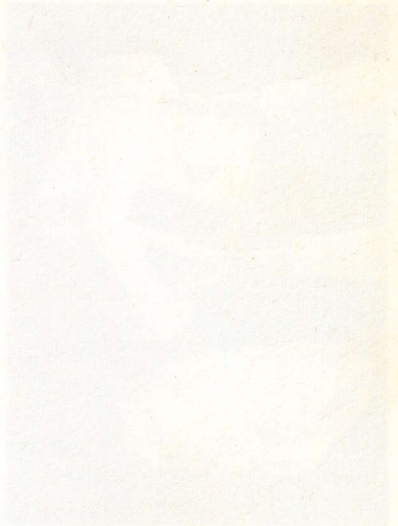
### Installation

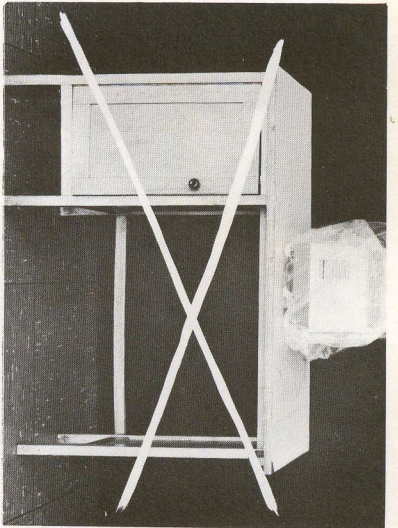Place balance on a sturdy table.  
Remove dust cover.

3

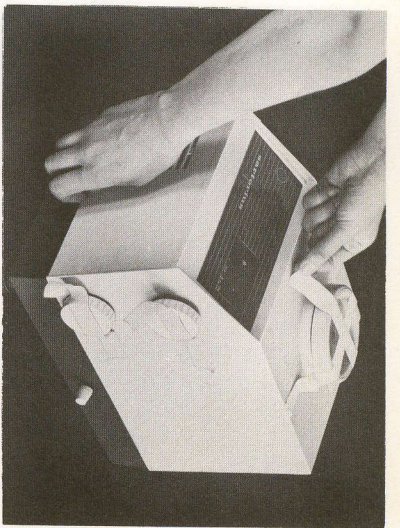Take off tape and remove front panel.

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
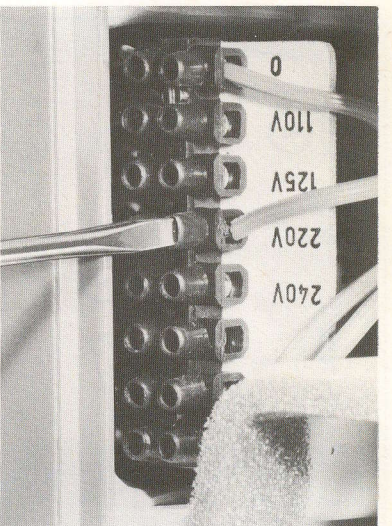
5

Make sure that balance is set for correct operating voltage. If necessary, change setting before connecting balance to line.

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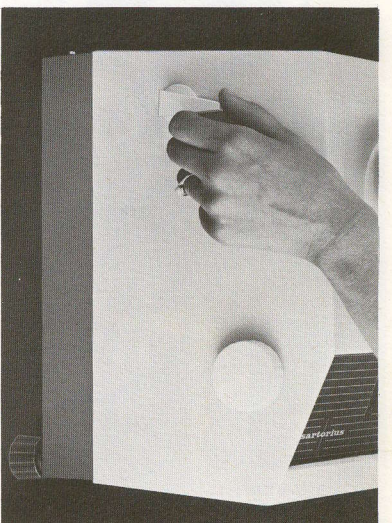
6

Place pan on support.  
Release transport arrestment.

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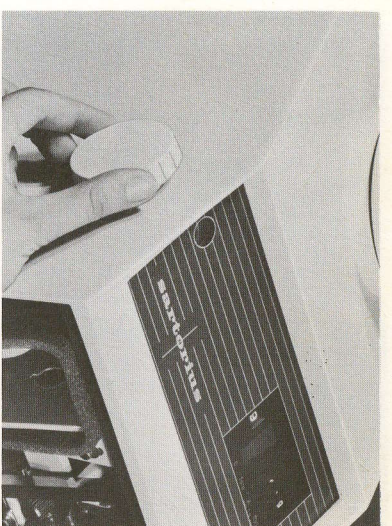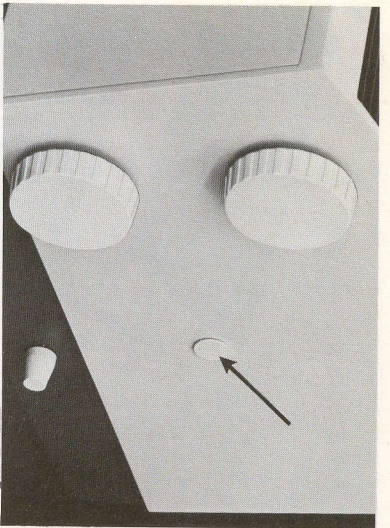
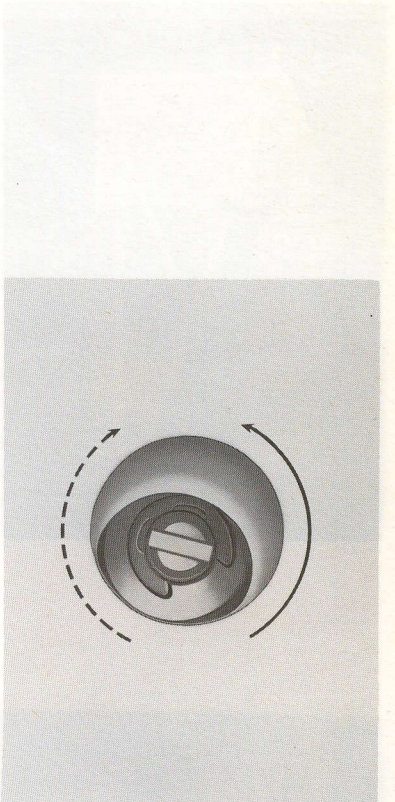
Tirer le bouchon.

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9

Loosen screw of transport arrestment with attached screw-driver.

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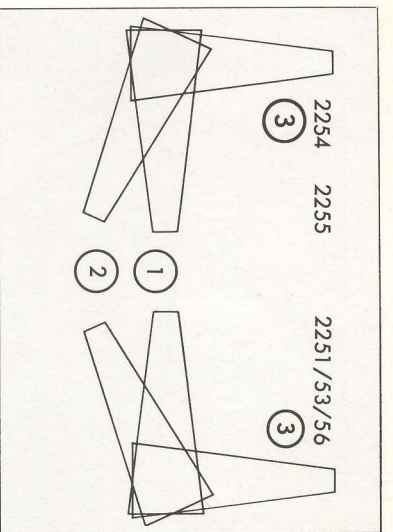
For transport, please fix screw again.

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
10

Balance in operation:  
 Balance shut off:  
 For transport only:

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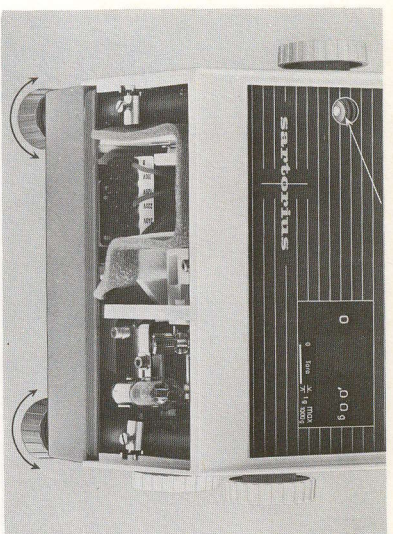
11

Level balance by means of foot screws and spirit level.

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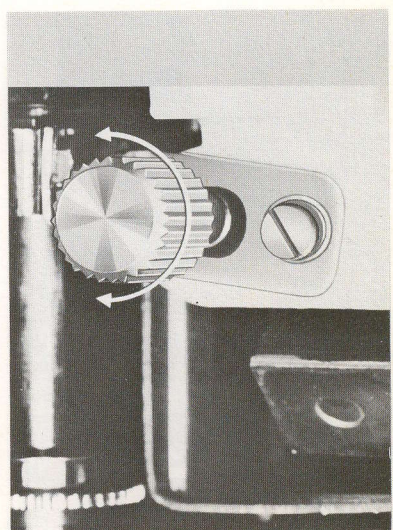
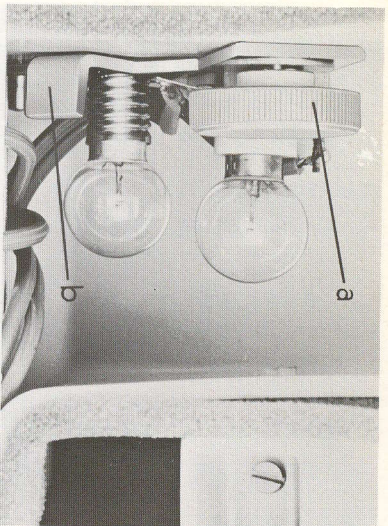
13

Adjust light intensity: Loosen knurled nut (a) re-position lamp bracket (b), re-tighten nut and re-insert front panel.

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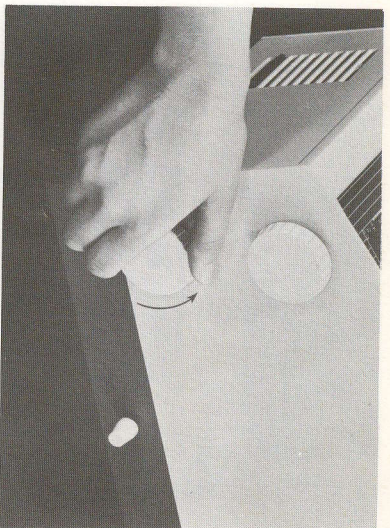
14

Turn tare knob to stop, then go back approx. 1/2 scale division.

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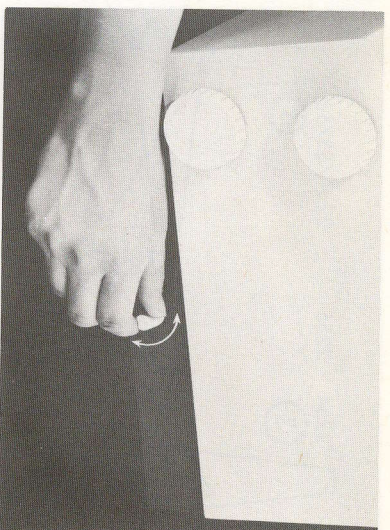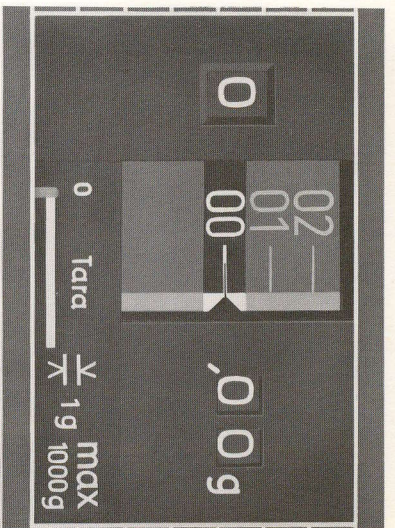
Ajústese el punto "0".

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
16

Zero point adjustment:  
Weight knob, micrometer and tare indicator  
must be at "0".

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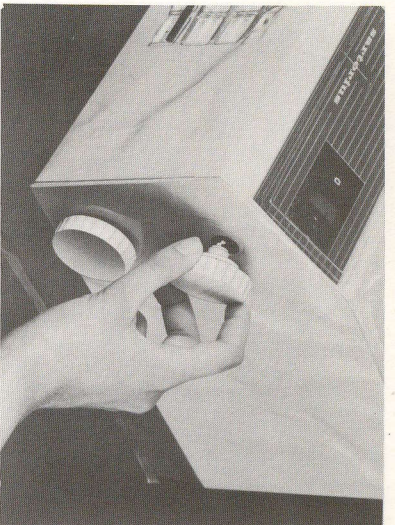
17

Pull off all operating knobs and levers, slip  
protective cover over housing and put knobs  
and levers back in place.

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
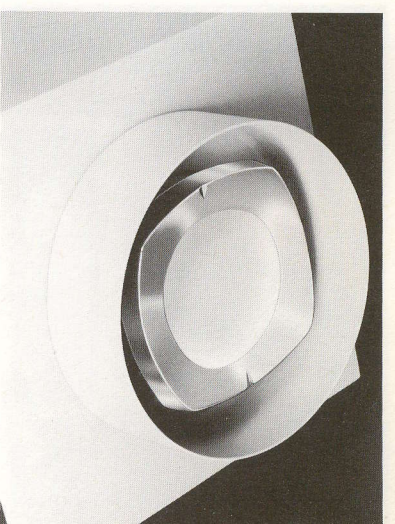
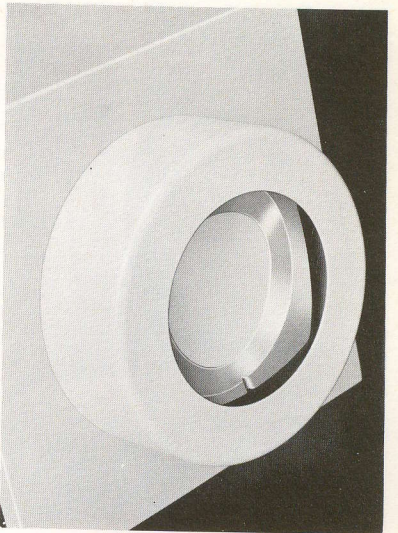
19

Model 2255 : draft shield is turned around  
for small beakers or samples.

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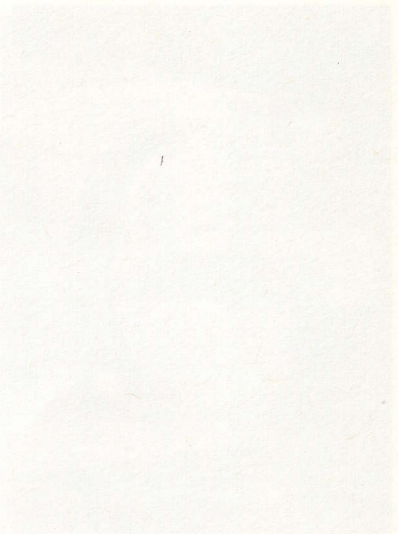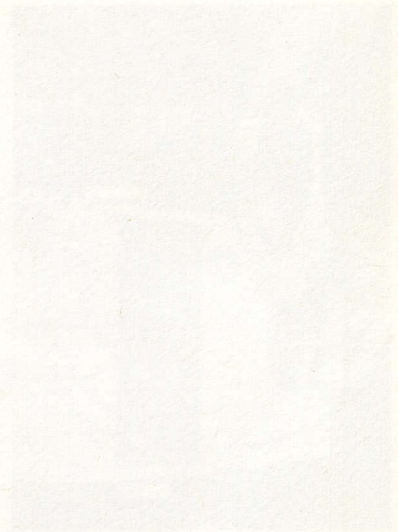
**Checking**

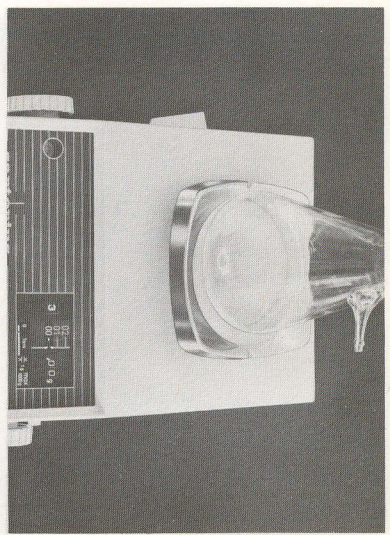Even though every balance leaves the factory in perfectly adjusted condition, slight changes may occur in transit.

20

Place an object of greater weight than the optical scale range on the balance. Set weight and tare knobs to "0".

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
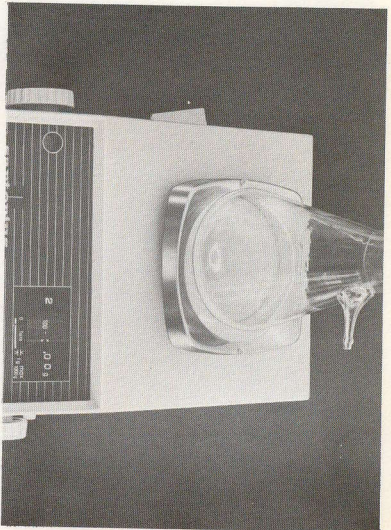
21

Return weight knob by 1 digit, scale must "swing in" on correct value.

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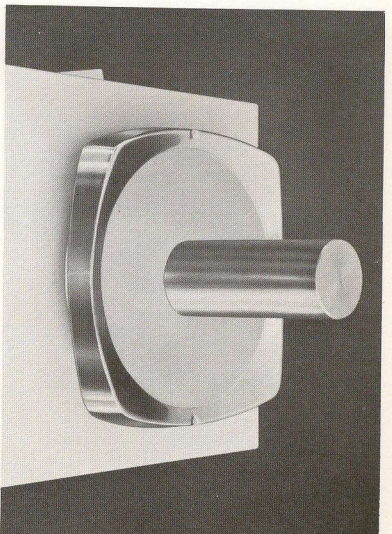
22

Model 2256 can be checked with 1 kg weight only (available on request).

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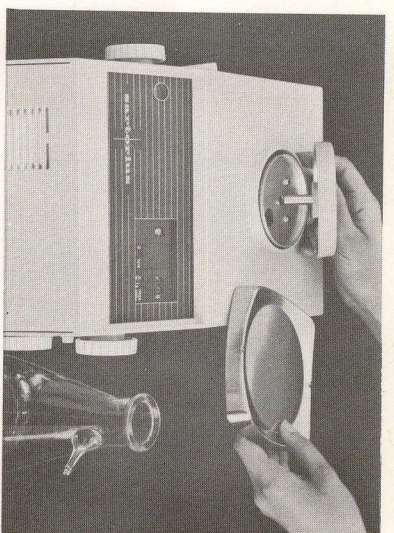
23

Remove pan and pan support if indication is incorrect.

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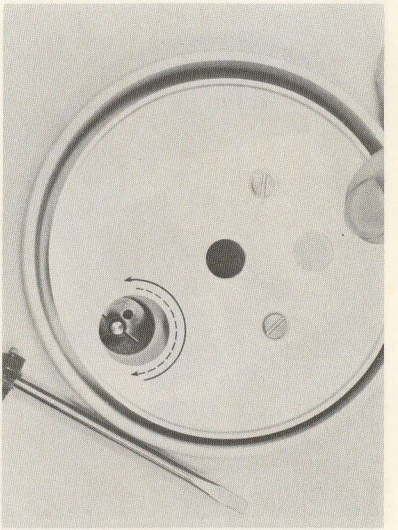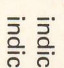
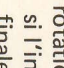
24

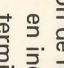
Models 2254/55: Remove stopper from balance housing. Push plate down.



Direction of rotation of the correcting weight :  
 indication below maximum scale deflection  
 indication above maximum scale deflection.

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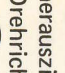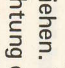
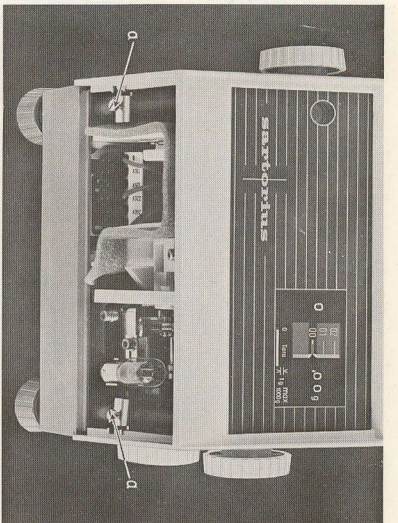
25

Models 2251/53/56 : Loosen screws (fig. a) and clamps on the rear panel. Pull off knobs and take off the casing.

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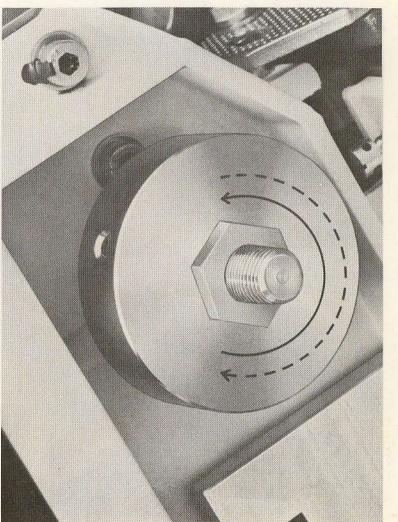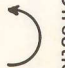
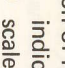
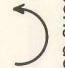
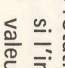
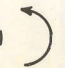
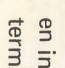
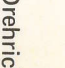
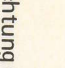
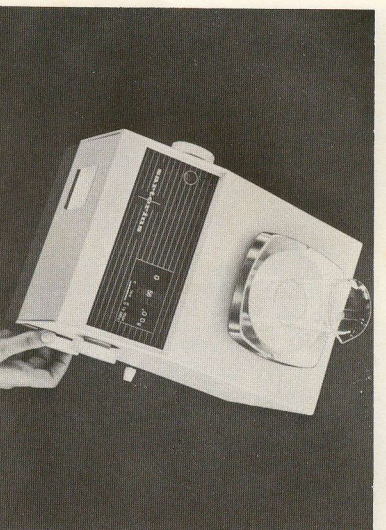
27

For taring, any value within the scale range can be tared to zero.

### Taring

The motion of the scale can easily be observed. No special indicator necessary.

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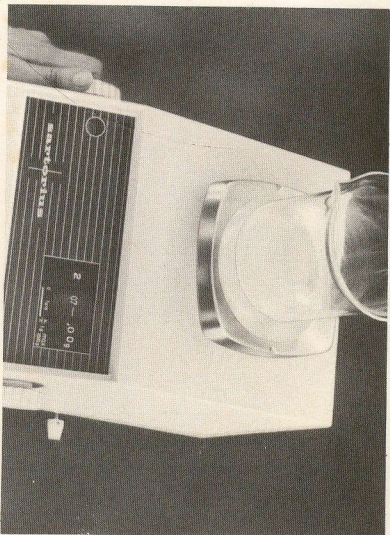
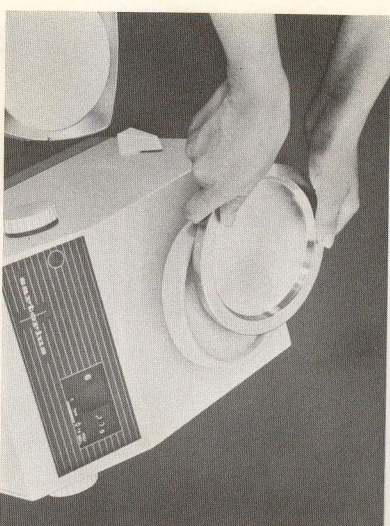
29

The tare range can be further extended if necessary by taking off the pan.

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
30

Model 2256: An additional approx. 2 kg tare weight t can be removed from below the pan.

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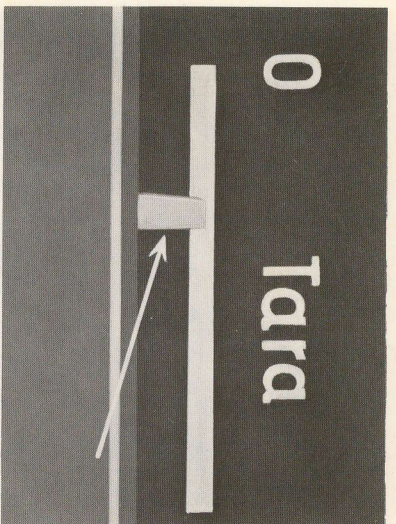
31

Pointer indicates taring position.

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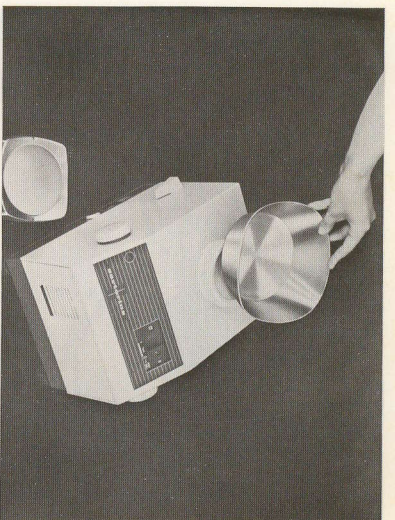32

Various size bowls (as special accessory) of matched weight are available (except for Model 2251).

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**Weighing to pre-set value**

Scale motion can be observed without special indicator.

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
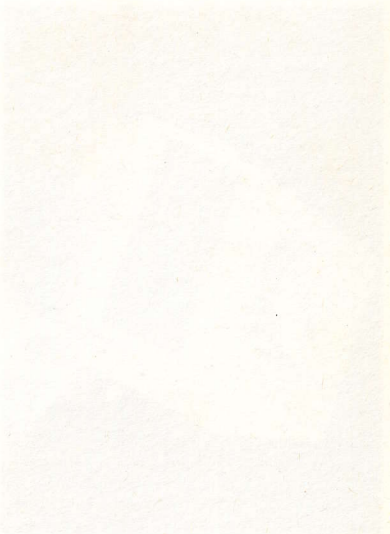
33

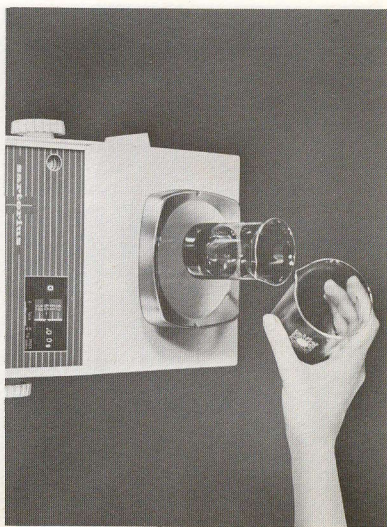
**Example: Sample weight 60 g.**  
After applying tare to compensate for weight of container, allow scale to swing to "60 g".

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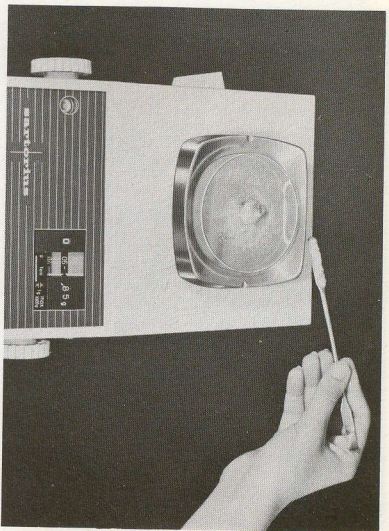34

**Example:** Sample weight 6.85 g (2254).  
Apply tare to compensate for container weight.  
Set micrometer to "85", allow scale to swing  
to "06".

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

### Weighing, regular

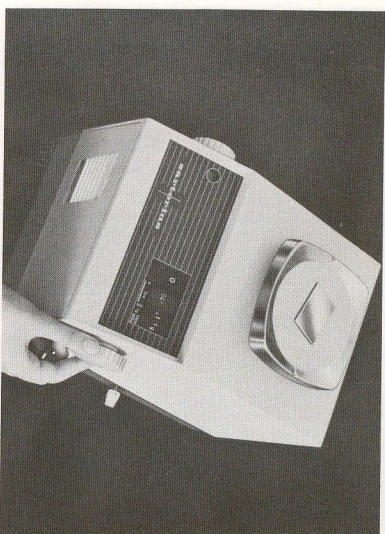35

Check zero adjustment. Place a load on the pan. Set micrometer after scale stops oscillating and read result.

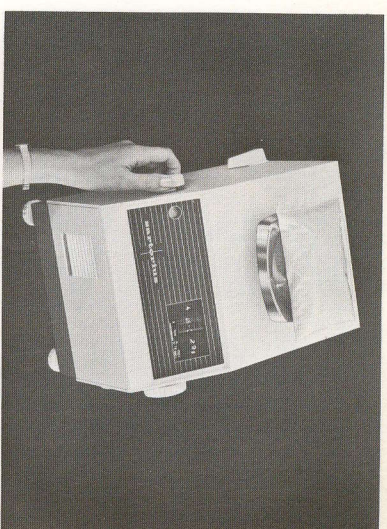36

If the weight of the sample is in excess of the optical scale range, turn weight knob until the scale "swings in", adjust micrometer and read result after scale stops oscillating.

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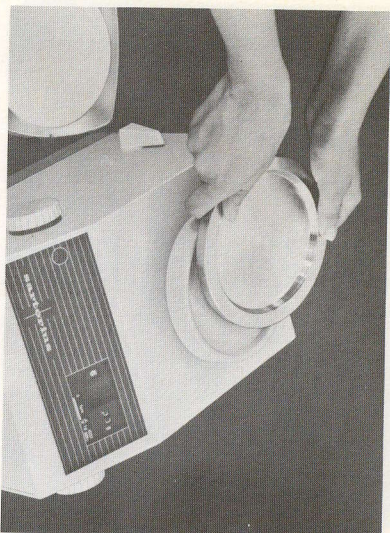
37


**Model 2251:** After removing the 2 kg-weight below the weighing pan, 2 kg are to be added to the weight-indication.

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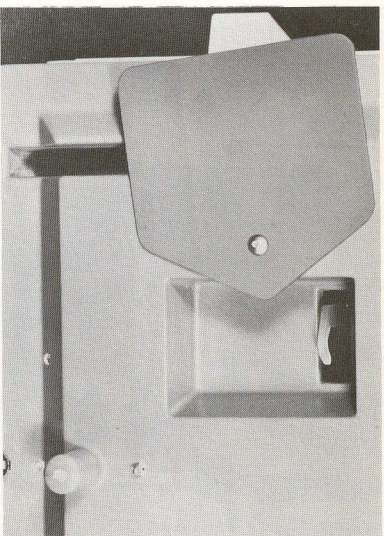38

### Weighing below the balance

All models are equipped with a hook for weighing below the balance. If necessary dial mechanical weights to maximum. Move lever (fig.6) into vertical position, place balance on one side and open cover sidewise.

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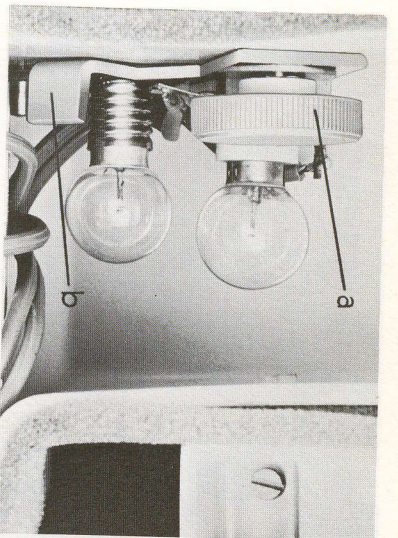
39

### Replacing the scale illuminator

Detach defective bulb and adjust light intensity according to fig. 13.

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### Zero point adjustment

Move tare knob (fig. 14) and zero adjustment knob (fig. 15) to halfway position. Unscrew the cover of the taring chamber (fig. 24) and add or remove taring material until the balance is at zero.

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**Preparing balance for transport**

Turn weight knob to maximum position,  
move lever (fig.6) up to vertical position.

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